


WOLFViewer 2.4 Download Instructions

If your WOLF Computer is Connected to the Internet:

1. If your WOLF computer is connected to the internet, you will receive automatic notifications to update WOLFViewer. This new built-in ability from 2.2 will ease your access to new software, avoiding on-site visits or remote connections with NanoCollect's Support team.
2. The next time you log into WOLFViewer, the software will prompt you to update. Click "yes" to continue.
3. WOLF2.4 will automatically be downloaded.
4. You should now see the new version in the WOLFViewer directory. Once you do, right click on WOLF2.2.200 and select delete.
5. WOLFViewer will close when the download is complete, and you can then launch the new version and are free to continue the cartridge set-up process as you normally would.

Example notification:

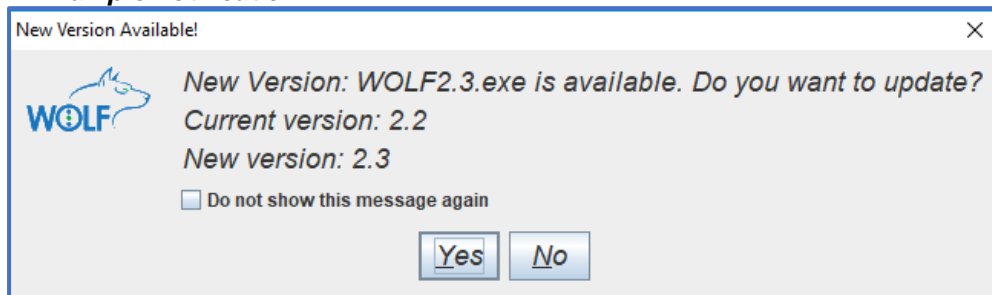

If your WOLF Computer is NOT Connected to the Internet:

1. Login to the [WOLF Users Knowledge Base](#).
2. Go to WOLFViewer Software section and click on the link to "WOLFViewer Software Version 2.4."
3. Download the file to a flash drive or another location to transport to your WOLF computer.
4. On the WOLF computer, paste the 2.4 .exe file into the WOLFViewer shortcut folder on your desktop where the current 2.2.200 version resides.
5. Right click on WOLF2.2.200 and select delete.
6. Utilize WOLF2.4 for all experiments moving forward.

Please contact your local FAS or Technical Support with any questions or issues.
support@nanocellect.com or 877-745-7678 (ext 3)